# NT-NK 4/1 AC

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# NT-NK 4/1 AC

## Power Supply for Low-Pressure- Mercury-Lamp NK 4/1

### 1. PACKING LIST

Power Supply  
Power Cord  
Lamp Cord

### 2. SPECIFICATIONS

<b>Lamp Current:</b>	~ 20 mA AC
<b>Lamp Voltage:</b>	~ 200V AC
<b>Ignition Voltage:</b>	>1000 V
<b>Power Input :</b>	230 V AC; 50 Hz
<b>Fuse:</b>	T200 mA
<b>Dimensions:</b>	W 11cm H 15cm D 35cm
<b>Weight:</b>	ca 3 kg

### 3. DESCRIPTION

The power supply is designed for the use of Heraeus Mercury lamp NK 4/1 in AC mode with a lamp current of about 20 mA. The usage of other arc lamps is also possible if the technical requirements meet the specifications of the power supply.

### 4. INSTALLATION

The lamp has to be connected to the supply by qualified personnel only.

**Attention : Since there is high voltage and the ignition voltage is over 1000 volts ensure enough insulation or space against physical contact to the lamps' electrodes !**

The attached lamp cord is for fixed installation into a lamp housing. Please regard for pin assignment to next page.

Ensure that the voltage value given on the back panel matches your power input and that power switch is off. Connect the power supply with the attached power cord to the socket. After switching power on lamp will ignite automatically.

## 5. PIN ASSIGNMENT

The lamp connector (3+1-poles) has this pin assignment

Pin	Assignment
1	sensor
2	lamp electrode
3	lamp electrode
	protective earth

Pin 1 is for detection if lamp is connected. For this ensure that pin 1 and 2 are only connected together if a lamp is installed.

Without a closure of pin 1 and pin 2 the usage of the lamp is impossible

**Caution! High Voltage at lamp socket!**

Service on lamp socket is permitted by qualified personnel only !

Disconnect the device from power before servicing the lamp socket !

## 6. SECURITY ADVICE

To prevent fire or shock, do not expose the unit to rain or moisture. The device should not be exposed to direct sun radiation. To avoid electrical shock, do not open the compartment.

Maintenance and repair service is permitted by qualified personnel only !